

### Additional Items Included But Not Shown:

PN 10337 Hex Key 2mm (for pot knob) PN 10490 Hex Key 5/64 inch (for brass tip set screw) PN 10566 Hex Key 2.5mm (for 3mm screws) PN 10610 Cable Tie 6in (3pcs)



В

K-

А

PN 10733 Power Supply - XP Power 5V 1A and PN 10734 Cable Assy 1.35mm ID x 3.5mm OD RA plug to USB A, 6 foot.



Tape to secure wires.



Power supplies are substitution without availability issues and changes in

PN 10736 Rubber plug to block unused AC power receptacle. Wayne Bonin 10/20/2021 PN 10976 Nikon Eclipse TS100 REV Installation Instructions 6 SHEET 1 OF 8

4

В

 $\rightarrow$ 

Nanodyne Replacement Illuminator for Nikon TS100 Microscope Installation Instructions: Step 1. Remove Illuminator/Stage Assy and Old Lamp Parts

3



(1) Remove the lamp assembly and stage from the microscope (three screws hold it in place).





(3) Remove the lamp holder assembly (three screws hold it in place).

This part will not be used.



(4) Remove the cover from the neck between the lamp housing and stage.

Remove the black plastic wire retainer (circled above) and the wires. Don't worry about damaging the retainer, it will not be used.

(2) Remove the black plastic housing by pressing firmly on two sides and lifting.

© Copyright 2021 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592

3

Δ

2



Wayne Bonin	10	/20/2	021
PN 10976 Nikon Eclips Installation Instructions	e TS1	100	REV 6
SHE	ΞТ 2	OF	8

Α

Nanodyne Replacement Illuminator for Nikon TS100 Microscope Installation Instructions: Step 2. Remove Pot Knob and Bottom Cover

3



4

(1) Remove the rubber cover from the intensity control knob to access the setscrew. This will take some force.



(2) Remove the metal knob by loosening the setscrew using the hex key provided.

Δ



(3) Remove the 6 screws circled above to remove the bottom cover.

Be aware that the the potentiometer shaft needs to be manuvered through the hole in the side of the base while removing the bottom cover.

© Copyright 2021 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592

3

 $\rightarrow$ 

Α

В

 $\left\{ + \right\}$ 

А

3 Nanodyne Replacement Illuminator for Nikon TS100 Microscope Installation Instructions: Step 3. Replace OEM Pot, Remove Old Wires



(1) Bottom cover removed - inside view.



(3) Install Nanodyne pot.



2

(4) Remove old wiring and route new pot cable, securing it as shown with wraps. Remove all connections to the AC power. Leave the green ground place.

© Copyright 2021 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-3

4

В

 $\rightarrow$ 

Δ

2

(2) Remove OEM Pot.

h 3 tie I wire in	in Wayne Bonin 10/20			021
	PN 10976 Nikon Eclipse TS100 Installation Instructions		REV 6	
-323-8592	SHEET	4	OF	8
	4			

1

1

В

А

# Nanodyne Replacement Illuminator for Nikon TS100 Microscope Installation Instructions: Step 4. Route Pot Cable, Connect Adapter and Install Illuminator



(1) Feed the pot cable end with the small connector through the back of the neck ....

### NOTE - Photos 3,4 and 5 were taken after the microscope was returned.

The adapter is installed over the microscope lens assembly shown in (2).



(2) ... and up and out at the lens assembly.





Δ



(5) Connect the illuminator to the adapter with the four 2-56 screws. Pull any excess cable back into the microscope neck.

© Copyright 2021 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592

2



В

(3) Place the adapter over the lens assembly (shown in (2)), with the pot cable going through the hole as shown above. Fasten the adapter in place with the M3 screws.



\*Note: Revised illuminator slides into the adapter and use set screw to tighten it instead of bolted as shown in (5).



Wayne Bonin 10/20/2021 PN 10976 Nikon Eclipse TS100 REV Installation Instructions 6 SHEET 5 OF 8

В

Nanodyne Replacement Illuminator for Nikon TS100 Microscope Installation Instructions: Pot Cable Connection Details

Connect the plug at the end of the Pot Cable Assembly to the mating socket of the illuminator, as shown in the pictures on this page. NOTE THAT THE PLUG IS KEYED TO ONLY GO INTO THE SOCKET ONE WAY, AS SHOWN.

Partially insert the plug into the mating socket of the illuminator by holding the wire next to the plug with your finger (photo 2).

Use your fingernails, if you have them, or tools like a tiny screwdriver or tweezers pushing on the side of the plug to fully insert it (photo 3).

The socket cannot be fully engaged by pushing on the wires, as the wires would just collapse.

To disconnect it if needed, pull the wire straight out by firmly gripping the black heat shrink tubing.







© Copyright 2021 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592

Δ

## Nanodyne Replacement Illuminator for Nikon TS100 Microscope Installation Instructions: Step 5. Complete Routing of Pot Cables



(1) Replace the bottom cover. Route the potentiometer cable through the slot in the cover as shown above.



(2) Re-attach the stage/neck/illuminator assembly to the microscope, and route the potentiometer cable up the back side of the scope and insert the loose end into the hole where the wire retainer used to be.

Use the VHB tape provided to secure the pot cable to the back of the microscope.

Connect the two pot cables as shown at left.



(3) Replace the cover on the neck as shown above.

Δ

2

© Copyright 2021 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592 3

4

4

В

 $\rightarrow$ 

А

Wayne Bonin	10/20/2021		
PN 10976 Nikon Eclipse Installation Instructions	TS1	00	REV 6
SHEET	7	OF	8

В

 $\left\{ +\right\}$ 

Α

Nanodyne Replacement Illuminator for Nikon TS100 Microscope Installation Instructions: Step 6. Replace Pot Knob, Install AC Power Blocking Plug.

3



4

В

 $\rightarrow$ 

(1) Replace the potentiometer knob and secure it with the hex key provided.

Replace the rubber covering, making sure to align the protrusion on the inside of the cover with the hole in the metal knob.



(2) Install PN 10736 rubber blocking plug in the receptacle on the back of the microscope as shown above (photo from generic microscope).

(3) Connect the USB end of the power cable to the power supply. Connect the other end of the power cable to the illuminator. Plug the power supply into an AC outlet and the illuminator is ready for use.

© Copyright 2021 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592 3

4

2

Wayne Bonin	10/20/2021		
PN 10976 Nikon Eclipse TS100 Installation Instructions		REV 6	
SHEE	т 8	OF	8

В

 $\leftarrow$ 

Α