


Nanodyne Replacement Illuminator for Olympus BX51-61 Bottom Illuminator: Included Items

Additional Items Included But Not Shown:

PN 10456 Hex Key 1.5mm (for microscope knob)
 PN 10697 Cable Ties.

PN 10407 BX50-51 Bottom illuminator Assy

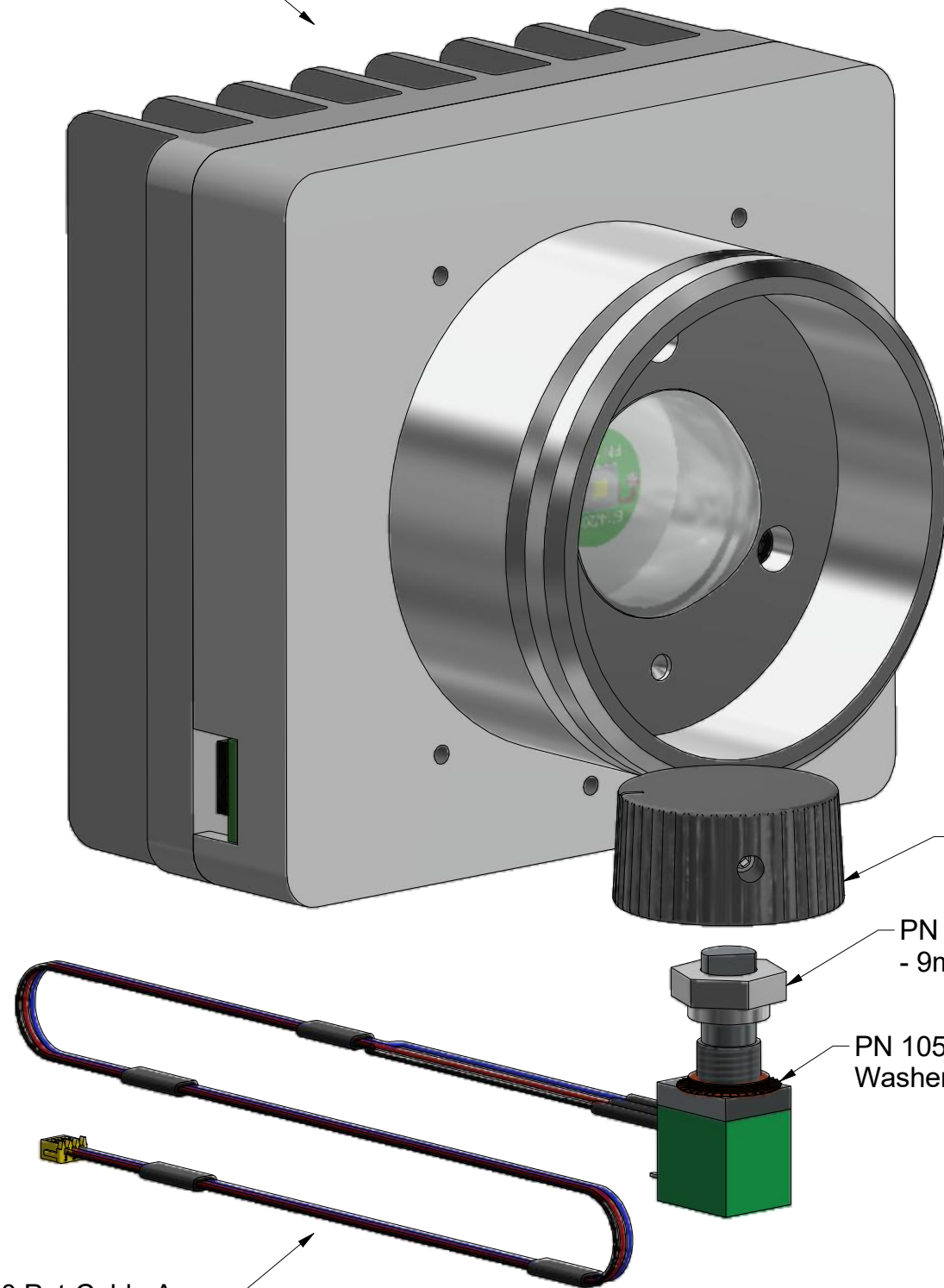

Pot Knob from BX51/61
(save from original microscope)

PN 11185 Custom Nut M7 x 0.75mm
- 9mm Bushing - 0.4375 in Hex

PN 10578 Serrated
Washer Assy

PN 11170 Pot-Cable Assy
(detabbed & shaft cut 5mm, 15 inch long)

The illuminator may be powered by plugging the cable into the power supply provided, or into a suitable USB port on a computer or other device.



3M VHB tape to secure pot cable



PN 11163 Power Supply 5V 2.1A and
 PN 10734 Power Cable 1.35mm ID
 3.5mm OD x USB A 6 ft.



PN 10736 Rubber plug to block unused
 AC power receptacle.

Power supplies are subject to change
 due to availability and regulations.

Wayne Bonin	7/8/2022	
PN 10413 Olympus BX51-61 Bottom illuminator System		REV 3
	SHEET 1	OF 1