

Nanodyne Replacement Illuminator for Nikon THP Pointer Light Installation Instructions: Step 1. Remove Old Bulb and Covers





Remove the bulb cover. It twists off by hand. The screwdriver is for later.



Remove the bulb.



В

Bulb after removal.





Remove the 4 screws holding the black rectangular cover using a phillips screwdiver. Remove the cover.

The screws and cover will be re-used.

Wayne Bonin	3/10	6/20	19
PN 11102 Nikon THP Pointer Light Installation Instruction			
	SHEET 2	OF	5


© Copyright 2019 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592

3

1

Nanodyne Replacement Illuminator for Nikon THP Pointer Light Installation Instructions: Step 2. Remove Obsolete Components



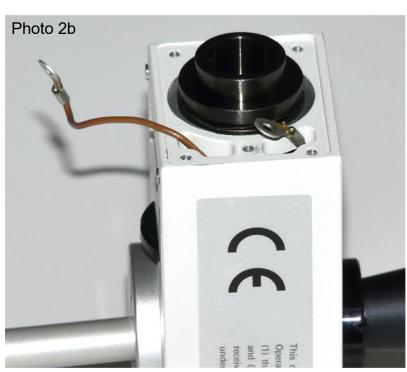

Remove the 2 screws holding the red wire to the sheet metal contact and the brown wire attached to the black cover



Re-attach the red wire terminal as shown in photo 2c to keep it out of the way. The brass contact will not be used. Cut the brown wire. It should be insulated with heat shrink tubing or electrical tape (not shown).





Remove the screw shown above to remove the brass lamp holder.

	P
and the second s	Photo 2e

Brass lamp holder removed from black cover. The lamp holder and screw are not used.

Wayne Bonin	3/16/20	19
PN 11102 Nikon Light Installation		REV 2
	SHEET 3 OF	5

© Copyright 2019 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592

Nanodyne Replacement Illuminator for Nikon THP Pointer Light Installation Instructions: Step 3. Attach Nanodyne light to OEM Cover



Nanodyne pointer light. Note the 2 tapped holes (one is partly hidden).



Note correct orientation of pointer light to cover.



provided. Tighten them with the 1.5mm hex key provided.





Ready to attach the cover with the Nanodyne pointer to the THP body.



Nanodyne pointer light installed.



Pointer light attached to cover.

Note that the illuminator is not centered on the cover, and the cover only fits one way on the THP body.

В

Be sure to attach the Nanodyne illuminator to the cover as shown in photo 3d, so the adjustment knob is in the front with the other controls (see next page).

> Wayne Bonin 3/16/2019 PN 11102 Nikon THP Pointer REV **Light Installation Instruction**


> > SHEET 4 OF 5

© Copyright 2019 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592


Nanodyne Replacement Illuminator for Nikon THP Pointer Light Installation Instructions: Step 4. Block Old Power Receptacle





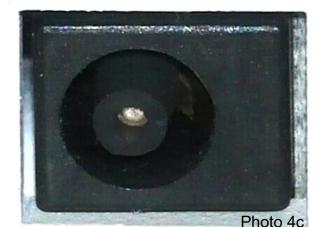

Closer views of blocked old receptacle.

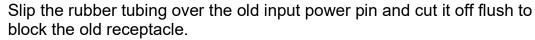



The power is applied by plugging in the new Nanodyne power cord to the receptacle shown in photo 4d between the white text.



Completed unit ready to attach to microscope.

Wayne Bonin	3/16/2019		
PN 11102 Nikon Light Installation	THP Pointer Instruction	REV 2	
	SHEET 5 OF	5	

В

© Copyright 2019 Nanodyne Measurement Systems. Document authorized for installation of Nanodyne equipment only. www.nano-dyne.com Wayne Bonin 651-323-8592